



## OVERVIEW

Fluo-Break is an emulsion breaker product from Emulseo. Fluo-Break has been formulated to efficiently disrupt droplets stabilized with Emulseo's surfactants. Fluo-Break allows an easy recovery of the droplet phase and is adapted for use with Emulseo's surfactants and oils.

## BENEFITS



Easy-to-use



Efficient recovery of the dispersed phase

## HOW TO USE

For best results:

- Collect the emulsion in a micro-centrifuge tube
- Carefully remove the excess oil from the bottom continuous phase
- Add Fluo Break to the emulsion (1 time the emulsion volume)\*
- Manually shake the micro centrifuge tube and centrifuge for 1 minute
- Recover the aqueous phase



\*Increasing the volume might be necessary for droplets stabilized with high concentration of surfactant.

## SAFETY NOTES

Wear protective gloves/protective clothing/eye protection.

For additional information, see the Material Safety Data Sheet.

## STORAGE

The emulsion breaker from Emulseo can be stored at room temperature in the supplied vial for six months without deterioration in performance.

## CONTACT

If you have any queries, please do not hesitate to e-mail us at: [contact@emulseo.com](mailto:contact@emulseo.com)